


Optics11

21-09-2017

OP1550 Interferometer driver issue:

A recent Windows update has caused an issue with running of the Piuma Nanoindentation system. It can be most easily recognised by the system not being able to perform a wavelength scan from the Calibration menu in the PC software. If after pressing the button, highlighted in the picture below, you get no result on the LCD screen of the OP1550 interferometer, you should follow the steps indicated below to solve the issue:

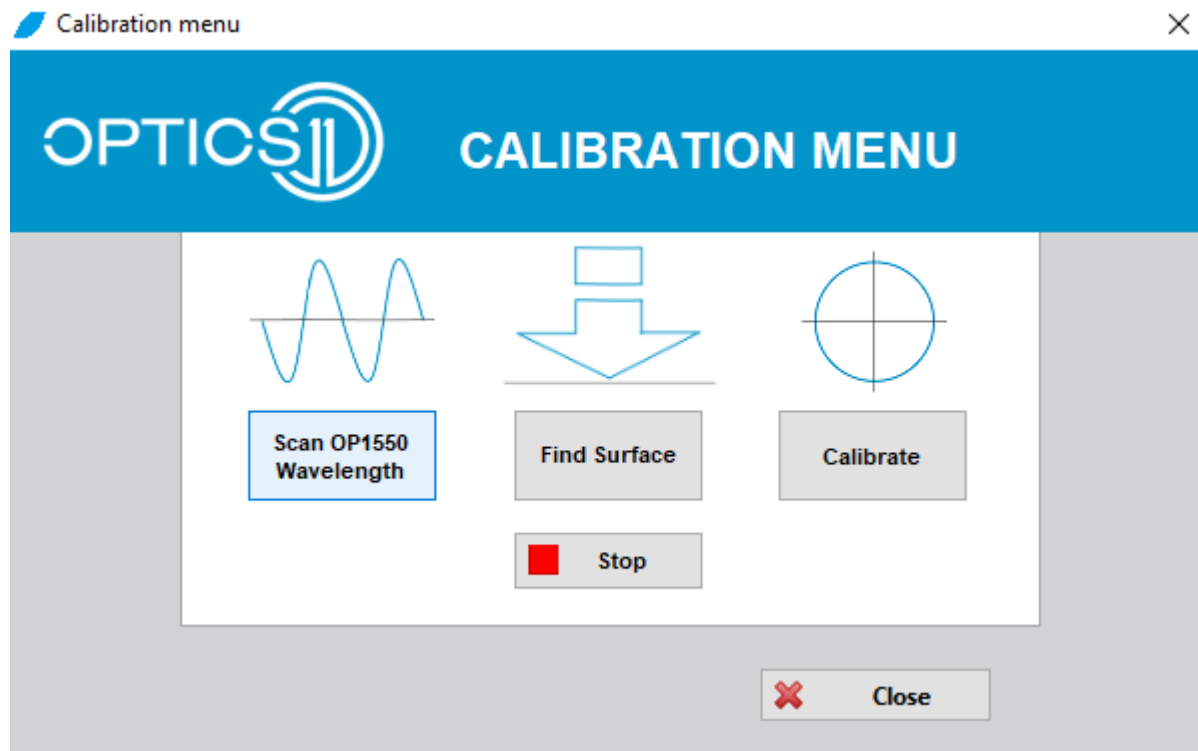

1. Download this file: [OP1550_Driver_Optics11.zip](#).
2. Extract the zip file, in the folder right click the 'Optics11InterferometerUsbDriver.inf' file. In the menu click 'Install' and go through the installation process.
3. The system should now be working again